## Ethidium homodimer 1, red fluorescent nucleic acid stain

<http://de.lumiprobe.com/p/ethidium-homodimer-1>

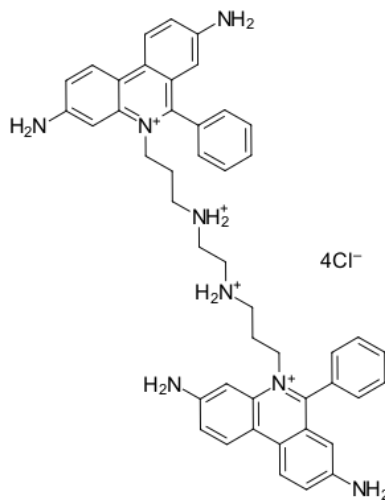
Ethidium Homodimer 1 (EthD-1) is a high-affinity, red fluorescent nucleic acid dye impermeant to live cells and widely used to identify dead cells in biological research.

EthD-1 is a dimeric analog of ethidium bromide, consisting of two ethidium moieties linked by a rigid bridge. Due to its large size and charged (dicationic) structure, EthD-1 cannot enter living cells with healthy membranes. In cells with compromised membrane integrity (dead or late-stage apoptotic cells), EthD-1 passively enters the cytoplasm and nucleus, where it intercalates into double-stranded nucleic acids (DNA and RNA) and exhibits a massive fluorescence enhancement (>40-fold) upon binding.


The EthD-1 red emission is easily distinguishable from green fluorescent probes (e.g., GFP, FITC, Calcein, [LUCS® 9](#), [LUCS® 13](#)), blue [Hoechst](#) stains, and far-red dyes ([LDS 751](#), [LUCS® 5](#)), enabling easy multiplexing.

EthD-1 is a standard dead-cell indicator in the viability and cytotoxicity assays, most commonly paired with green fluorescent live-cell stain like Calcein AM, [LUCS® 9](#), or [LUCS® 13](#). It is also used as a counterstain in fluorescence microscopy.

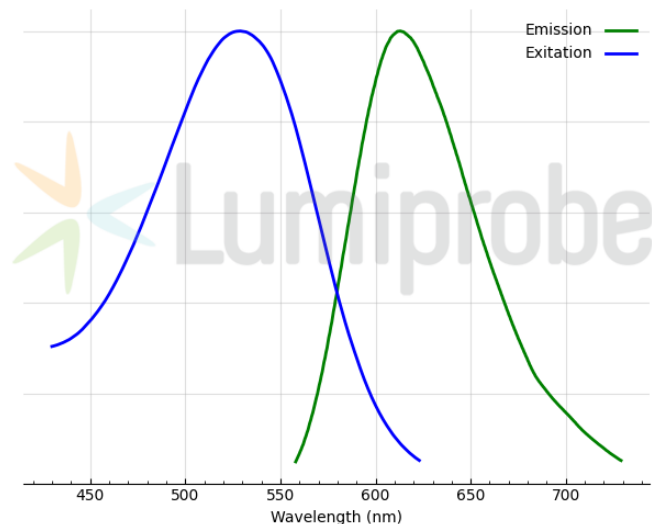Note that, while highly specific for dead cells in short-term assays, prolonged exposure can lead to its uptake by live cells, as it is mildly cytotoxic and can eventually permeabilize membranes.

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